pH-Fix 1.7-3.8





Please refer to instructions for use

REF 92190

Pack content:

100 test strips 1 reference color scale pH 1.7–3.8

Instructions for use:

- 1. Dip test strip for 1 s with all test fields in the sample solution.
- 2. Shake off excess sample solution.
- 3. Compare with color scale and read off the corresponding pH-value.



Note:

If sample solution is weakly buffered leave test strip in solution until color no longer changes.

Storage:

Avoid exposing the test strips to sunlight and moisture. Store the container in a dry place (storage temperature 4-30 °C).

Disposal:

Used test strips may be disposed of as domestic waste.



Please read instructions for use!



Item number



Manufacturer

en