

# pH-Fix 5.1–7.2

## Test strips for measuring pH values in the range of 5.1–7.2



Please refer to instructions for use

REF 92140

### Pack content:

100 test strips

1 reference color scale pH 5.1–7.2

### Instructions for use:

1. Dip test strip for 1 s with all test fields in the sample solution.
2. Shake off excess sample solution.
3. Compare with color scale and read off the corresponding pH-value.



### Note:

If sample solution is weakly buffered leave test strip in solution until color no longer changes.

### Storage:

Avoid exposing the test strips to sunlight and moisture. Store the container in a dry place (storage temperature 4–30 °C).

### Disposal:

Used test strips may be disposed of as domestic waste.



Please read instructions for use!



Item number



Manufacturer

