

# PEROXTESMO MI

en

## Test paper for the detection of lactoperoxidase in milk

The enzyme lactoperoxidase is always present in raw milk. It will only be deactivated by high temperature heating (UHT-milk) but not through normal pasteurization. So this test paper reacts with UHT-milk lactoperoxidase negative and with pasteurized milk lactoperoxidase positive.

### Presentation:

Box of 100 test papers 15 x 15 mm.

### Color reaction:

Peroxtesmo MI reacts with lactoperoxidase by a coloration of blue flecks on a white background.

### Limit of sensitivity:

3% raw milk in UHT-milk

### Instructions for use:





Remove one test paper from the box. Immediately close the box. Apply one drop of milk on the test paper and within 2 minutes the presence of lactoperoxidase will be detected by a blue fleck.

### Comment:

Any color change after 2 minutes should be ignored.  
The presence of Azidiol leads to a weaker color reaction.

### Storage conditions:

Protect from direct sunlight and moisture. Store container in a cool and dry place (storage temperature below 20 °C).

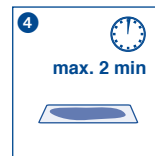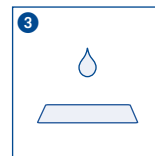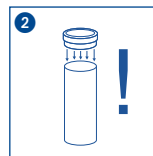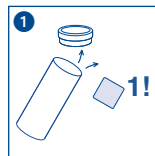

REF	Artikelnummer / Item number / Référence produit / Referencia / Artikelnummer / Numero di catalogo / Numer artykułu	LOT	Chargencode / Batch identification / Número de lot / Código de lote / Productienummer / Codice del lotto / Numer partii
!	Achtung! / Warning! / Attention! / ¡Atención! / Let op! / Attenzione! / Uwaga!		Verwendbar bis / Use by / À utiliser avant / Fecha de caducidad / Te gebruiken tot / Utilizzare entro (anno / mese) / Przydatność do użycia
	Temperaturbegrenzung / Permitted storage temperature range / Limites de température / Limites de temperatura / Temperaturbegrenzung / Limite de temperatura / Ograniczenie temperatury		Trocken aufbewahren / Store in a dry place / Conserver au sec / Mantener seco / Op een droge plaats bewaren / Mantener asciutto / Przechowywać w suchym miejscu
i	Gebrauchsanweisung beachten / Please read instructions for use / Respecter les instructions d'utilisation / Obsérvense las instrucciones de uso / Lees de bijsluiter / Consultare le istru- zioni per l'uso / Przestrzegać instrukcję użycia		Packung geschlossen halten / Keep container closed / Refermer la boîte / Mantenga el envase cerrado / Verpak- king gesloten houden / Conservare la confezione chiusa / Przechowywać pojemnik szczelnie zamknięty
Σ	Ausreichend für <n> Prüfungen / Contains sufficient for <n> tests / Contenu suffisant pour <n> tests / Contenido suficiente para <n> tests / Voldoende voor <n> tests / Contenuto sufficiente per <n> test / Wystarczająco dla kontroli <n>		